



**BIOPAC Systems, Inc.**

42 Aero Camino  
Goleta, CA 93117  
Ph (805) 685-0066  
Fax (805) 685-0067  
[www.biopac.com](http://www.biopac.com)  
[info@biopac.com](mailto:info@biopac.com)

---

## **iMac and G3 Compatibility Issues**

➡ See the [USB1M](#) Solution

---

This gPort<sup>®</sup> adapter solution was replaced with the USB1M

*BIOPAC Systems, Inc.* has a solution to compatibility issues for communication between the iMac and G3 line of Apple Computers and the MP30 and MP100 units.

Griffin Technology has introduced a number of compatibility products for the blue and white G3 Macs. The gPort<sup>®</sup> adapter provides standard Mac serial interface that is compatible with all Mac serial devices. It supports all Apple and third party serial and local talk printers, midi devices, as well as any other serial device that works on a Mac.

The gPort<sup>®</sup> uses the modem slot on your B&W G3 so it provides "board level" compatibility. It is also geoport compatible. The gPort<sup>®</sup> takes only minutes to install.



The iPort® universal adapter for iMacs has the I/O connections Mac users have come to depend on. Designed from the ground up with the iMac in mind, it features video out capable of supporting up to 1600x1200 resolution. It also supports video mirroring in iMac compatible video modes.

The iPort® has a serial port that is compatible with all serial devices that currently work with a Macintosh computer. It also supports local talk networking. You can connect to a local talk network and print or share files jut like Macs have always been able to do.

**The iPort® is only compatible with revision A & B "Bondi Blue" 233MHz and 266MHz iMacs.**

- The iPort® is not compatible with the 'C" series iMacs.
- Contact Apple Computer to find out which series iMac you own.



- Griffin Technology may have a compatible adapter for the new flavored machines. Please contact Griffin Technology for inquiries or questions:

Griffin Technology  
80 Fesslers Lane  
Nashville, TN 37210 USA  
Tel: 615 255 0990  
Fax: 615 255 8040  
E-Mail: [sales@griffintechnology.com](mailto:sales@griffintechnology.com)